

DuraUse™ Anti-Spoil Skim Milk Powder

3 lbs



15875 Gaither Drive
Gaithersburg, MD 20877
FAX. 301-560-4919
TEL. 301-330-5966
Email: info@signagenlabs.com
Web: www.signagenlabs.com

This product is for laboratory research ONLY and not for diagnostic use

Description:

This product is specially formulated for western blotting use; Nonfat milk powder (5 %) in TBST solution is utilized for blocking the unspecific binding in Western Blot. Incubate the membrane either PVDF or nitrocellulose with 5 % defat milk powder in TBST for 1 hour at room temperature or at 4 °C overnight.

Features:

This product is screened to guarantee good blocking effect with azide and antibiotics added; If prepared properly, able to store the prepared blocking solution (5 % milk powder in TBST) at 4 °C for 3 months.

Instruction:

1. Shake well before weighing.
2. Dissolve nonfat milk powder in TBST solution and make final concentration 5 %
3. Store the 5 % nonfat milk powder in TBST at 4 °C.
4. The nonfat milk powder in TBST can be preserved for 3 months with the sterilized container and appropriate preparation.

Storage: Upon arrival store this product at room temperature. For long term, store the product at 4 °C. Product shipped at ambient temperature.